

BioFX™ Tris Buffered Saline/Casein Block and Diluent-10X Concentrate

Product insert

Product number:

TBSC-0100-01 (100 mL)
 TBSC-1000-01 (1000 mL)
 Custom Sizes Available

Product use:

BioFX™ Tris Buffered Saline/Casein Block and Diluent-10X Concentrate minimizes non-specific interactions of the secondary reactants with each other and with the primary solid phase binding sites. The high blocking efficiency of casein is due to its content of small molecular weight proteins. The solution is ready for dilution for specific application requirements.

Product stability, storage and specifications:

Product stability	2 years from date of manufacture.
Storage	2-8°C
Specification	The solution should be clear and colorless to light yellow in appearance. To minimize the risk of contamination, a fresh working solution should be prepared each day or store unused working solutions at the recommended storage temperature.

Recommendations for use:

- 1) Dilute 1 mL TBS Casein Block/Diluent-10X Concentrate in 9 mL reagent quality water to produce a 1X working solution.
- 2) Allow the 1X working solution to equilibrate to room temperature (25°C) prior to use.
- 3) To use as a blocking reagent, completely cover the target surface with the 1X working solution. To use as a diluent, dilute the antigen/antibody to a desired concentration in the 1X working solution and completely cover the target surface as needed.
 - An overnight incubation step at 2-8°C following the coating process will yield the best results.

For technical assistance, email ivdtechsupport@surmodics.com

Related products:

In-Solution Protein Stabilizers & Diluents:
StabilZyme™ HRP Conjugate Stabilizer (SZ02)
StabilZyme™ SELECT Stabilizer (SZ03)
StabilZyme™ NOBLE Stabilizer (SZ04)
StabilZyme™ Protein Free Stabilizer (SZPF)
Surmodics™ Assay Diluent (Protein-Free) (SM01)
MatrixGuard™ Diluent (SM02)
Blockers/Stabilizers:
StabilGuard™ Immunoassay Stabilizer-BSA-Free (SG01)
StabilCoat™ Immunoassay Stabilizer (SC01)
StabilBlock™ Immunoassay Stabilizer (ST01)
BioFX™ Phosphate Buffered Saline/Casein Block and Diluent – 5X Concentrate (PBSC)

BioFX™ Tris Buffered Saline/Casein Block and Diluent-10X Concentrate

Product insert

Substrates:
BioFX™ TMB Super Sensitive One Component HRP Microwell Substrate (TMBS)
BioFX™ TMB One Component HRP Microwell Substrate (TMBW)
BioFX™ TMB Conductivity One Component HRP Microwell Substrate (TMBC)
BioFX™ TMB Slow Kinetic One Component HRP Microwell Substrate (TMSK)
BioFX™ TMB Extended Range One Component HRP Microwell Substrate (TMBX)
BioFX™ TMB Enhanced HRP Membrane Substrate (ESPM)
BioFX™ TMB One Component HRP Membrane Substrate (TMBM)
BioFX™ Chemiluminescent Ultra Sensitive HRP Membrane Substrate (CHMM)
BioFX™ AP Yellow One Component Microwell Substrate (pNPP) with Stabilizing Pellets (PNPS)

Products distributed by



DIARECT GmbH
Bötzing Str. 29 B
79111 Freiburg
Germany

Tel. +49 (0) 761 47979-0
Fax +49 (0) 761 47979 -29

www.bbisolutions.com
orders-dia@bbisolutions.com
info-dia@bbisolutions.com

SURMODICS, BIOFX, STABILZYME, MATRIXGUARD, STABILCOAT, STABILGUARD and STABILBLOCK are trademarks of Surmodics, Inc. and/or its affiliates.

These products are for further manufacture and/or to be used as a component with in-vitro diagnostics immunoassays and are not intended for use in human or therapeutics purposes. Sales are without any seller's warranty or representation, expressed or implied, by usage or otherwise; no claims beyond replacement of unacceptable material or refund of purchase price shall be allowed. All claims must be made according to DIARECT's Terms and Conditions of Sale.